

Quick Change Cuffs

Time is your labs most valuable commodity and Coy Labs is always looking for ways to minimize your anaerobic chamber maintenance. Now the latest innovation for the most cited anaerobic chamber in the world the **Quick Change Cuff (QCC)** system for the Coy Vinyl Anaerobic Chambers.

What use to be a 20-30 minute exchange is now a **30 second or less** with the QCC.

Exchange process is quick and easy.

1. A port plug with handle is used to seal the port while the old glove is removed.
2. A new glove attached to an adaptor ring is brought into the chamber.
3. Adaptor Ring is pushed on to the cuff port
4. Heavy rubber seals are rolled on to the port to secure the glove to the port cuff



Advantages

- Significant gas savings as damaged gloves are readily exchanged
- Appropriate size gloves are quickly exchange for any size user.



Ask your Coy representative about a gloveless version